

Mode

■ **list wave delay** :

Settings for list-pileup or list-wave mode.
Adjust the delay of the acquired waveform.
Set it near 5digit.
It can be set from 1digit to 31digit.

■ **list wave data** :


Parameters for list-pileup mode or list-wave.
Set the number of data points for pile-up waveform output.
Can be set from 200ns to 2μs in 8ns increments



Example of use

Model : APV8516-14

- ADC: 500MSPs, 16CH, resolution 14-bit
- Time resolution: Coarse 2ns | Fine 7.8ps, LSB
- Throughput: 1Mcps and more / ch.

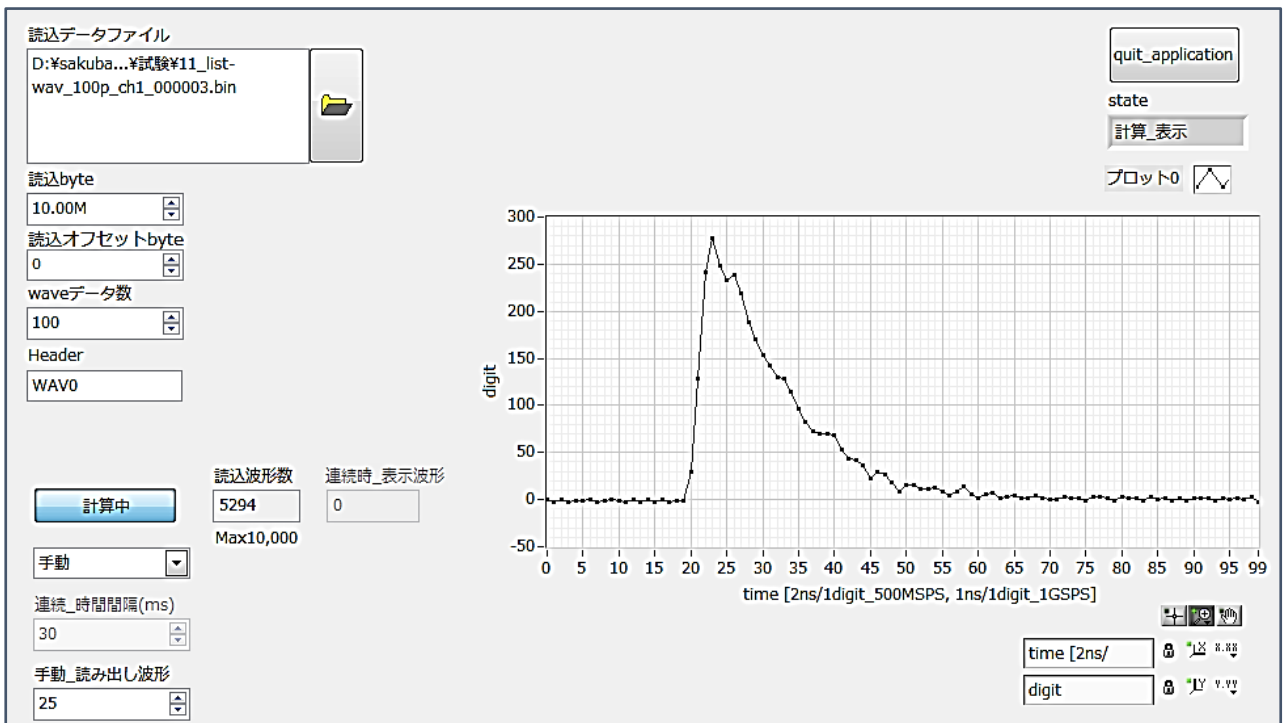

Waveform read by test application

*Images is for illustration purpose.

*Please note that contents may change without prior notice.

